

WaveTrain[®] 2

External Cavity Frequency Doubler For CW Lasers



WaveTrain 2 is an advanced stand-alone device that provides simple frequency doubling of single frequency CW laser beams. With greater efficiency than traditional intra-cavity or extra-cavity methods, this product can be used to generate the second harmonic of the output of the Matisse[®] 2.

WaveTrain 2 utilizes a ring cavity to achieve efficiencies greater than 35%. Unlike conventional bow-tie cavities, WaveTrain 2 is based on a patented triangle-shaped cavity configuration called the DeltaConcept[®], which allows the cavity length to be adjusted with no effect whatsoever on output beam position, direction or beam astigmatism.

Output stability is further enhanced by the use of a lightweight piezo-activated prism to optimize the cavity length, making the WaveTrain 2 especially resistant to vibrations and acoustic noise. This inherent beam stability enables the control hardware to automatically lock the cavity for maximum doubling efficiency during both fixed frequency and scanned operation.

Ease of access and the use of pre-mounted modular optics and crystals insure maximum wavelength flexibility, enabling the WaveTrain 2 to be operated effectively over the entire 410–1600 nm input wavelength range.

The WaveTrain 2 Advantage

- Input wavelength range 410 to 1600 nm
- Fast servo loop ensures power stability
- Continuous scans of 65 GHz at 10 GHz per second
- Sealed oscillator block with external adjustments prevent contamination of optics
- Only two pre-mounted mirrors for quick change of wavelength ranges
- Low cavity losses and high doubling efficiencies


Applications

- High-resolution UV spectroscopy
- Atom cooling and magneto-optic trapping
- Tunable UV experiments
- Bose-Einstein condensates
- Atomic clocks




WaveTrain 2


WaveTrain 2 Efficiency Performance^{1,2}

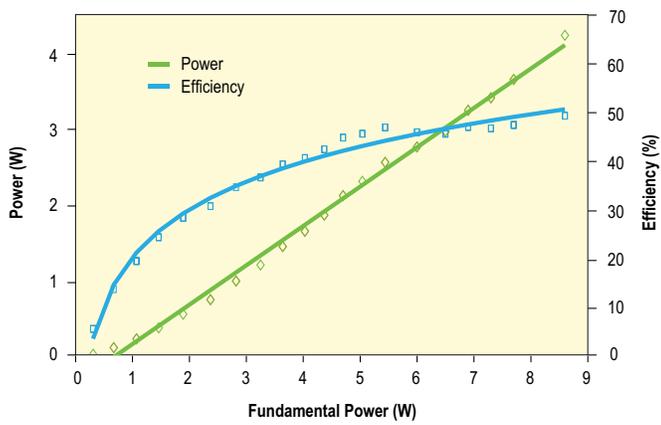

- Typically measured performance; not a guaranteed or warranted specification.
- Taken with LBO crystal at 754 nm.

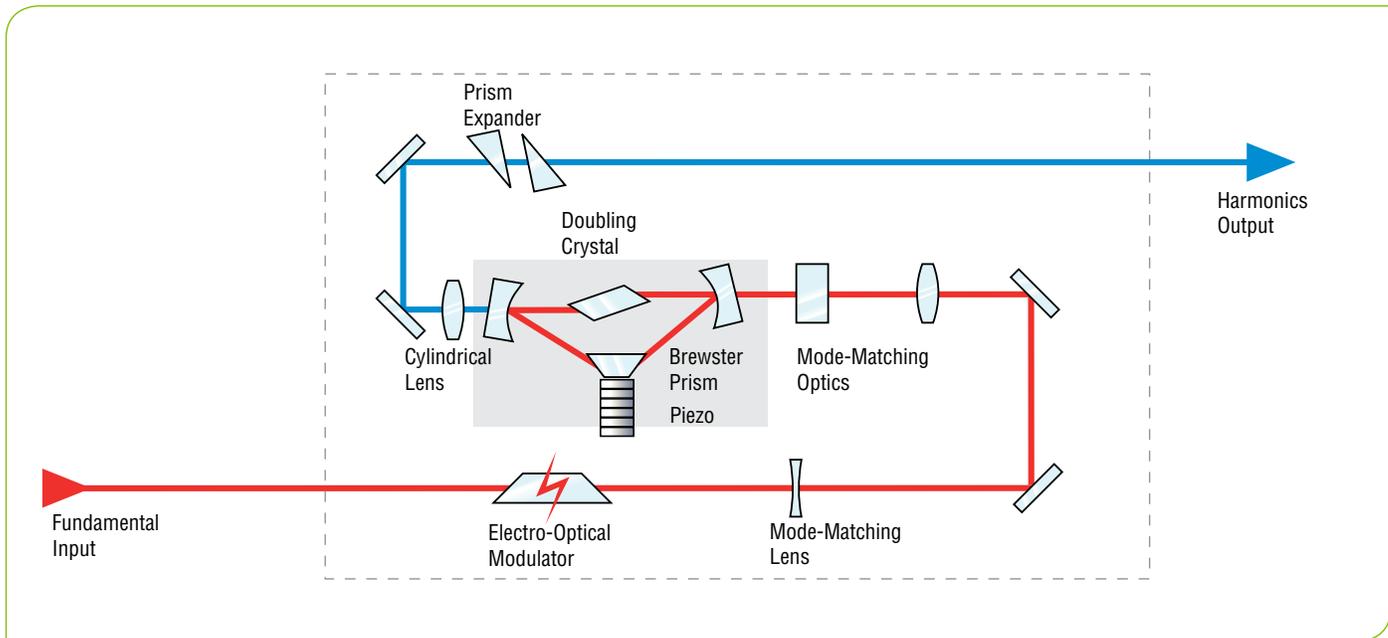

WaveTrain 2 Optical Layout

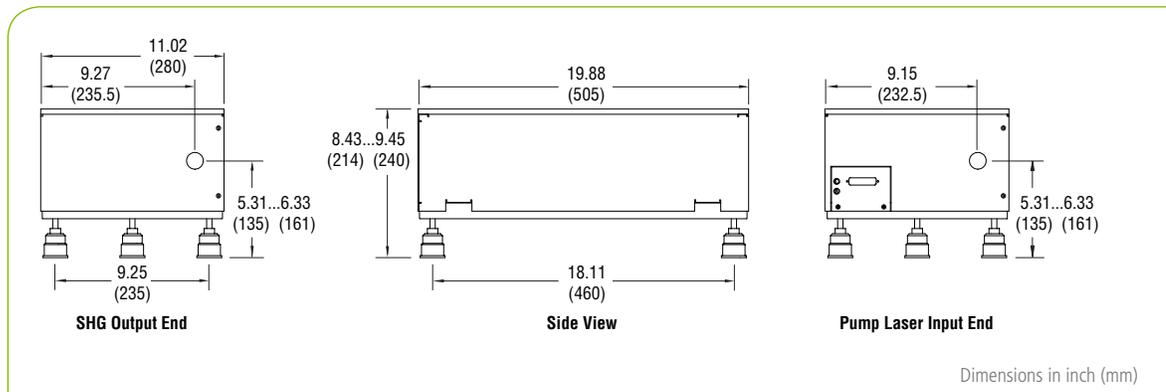

WaveTrain 2 Dimensions

WaveTrain 2 Specifications¹

Wavelength Range	670–1050 nm	550–780 nm	410–550 nm	1050–1600 nm
Automated Scan Range	>40 GHz	>65 GHz	>65 GHz	>40 GHz
Phase-Match Tuning ^{2,3}	±25 nm @ 800 nm	±15 nm @ 660 nm	Contact factory	Contact factory
Maximum Scan Rate	10 GHz/sec (fundamental)			

Efficiency⁴

Input Power	670–1050 nm	550–780 nm	410–550 nm	1050–1600 nm
300–600 mW	>6%	>4%	4–10%	>12%
600–1300 mW	>15%	>10%	4–10%	>12%
1300–3000 mW	>25%	>15%	4–10%	>12%
3000–10000 mW	>35%	>15%	4–10%	>12%

Requirements

Pump Laser	TEM ₀₀ , M ² <1.1, single frequency laser Linewidth <10 MHz, linear polarized
Ambient Conditions	Constant temperature in the 20–25°C range
Laboratory	Vibrational isolated optical table, dust-free air (flow box)
Voltage	115/230 V, single phase, 50/60 Hz

Physical Characteristics

Size (L × W × H)	19.88 x 11.02 x 8.43–9.45 in (505 x 280 x 214–240 mm)
------------------	---

1. Due to our continuous product improvement program, specifications are subject to change without notice.

2. Phase-match tuning is defined by an intensity reduction of 50% in the SHG.

3. Crystal cut wavelength may vary by 1%.

4. Specifications for 670–1050 nm and 550–780 nm tuning range only with Matisse Ti:Sapphire and dye lasers.



www.spectra-physics.com

1565 Barber Lane, Milpitas, CA 95035 USA

PHONE: 1-800-775-5273 1-408-980-4300 FAX: 1-408-980-6921 EMAIL: sales@spectra-physics.com

Belgium +32-(0)0800-11 257 belgium@newport.com

China +86-10-6267-0065 info@spectra-physics.com.cn

France +33-(0)1-60-91-68-68 france@newport.com

Germany / Austria / Switzerland +49-(0)6151-708-0 germany@newport.com

Japan +81-3-3556-2705 spectra-physics.jp@mksinst.com

Korea +82-31-8021-1600 korea@spectra-physics.com

Netherlands +31-(0)30 6592111 netherlands@newport.com

Singapore +65-6664-0040 sales.sg@newport.com

Taiwan +886-3-575-3040 sales@newport.com.tw

United Kingdom +44-1235-432-710 uk@newport.com